


LED BINOCULAR MICROSCOPE 60X

INSTRUCTIONS:

1. Place the sample to be viewed on a flat surface.
2. Place the microscope vertically over the sample with the Objective Lens (D) facing downwards.
3. If additional light is needed, turn on the LED lights using the Switch (F).
4. View the magnified object through the Eye Lens (A).
5. If the image appears out of focus, use the Focusing Barrel (C) to make adjustments.
6. Turn off the LED lights when finished.

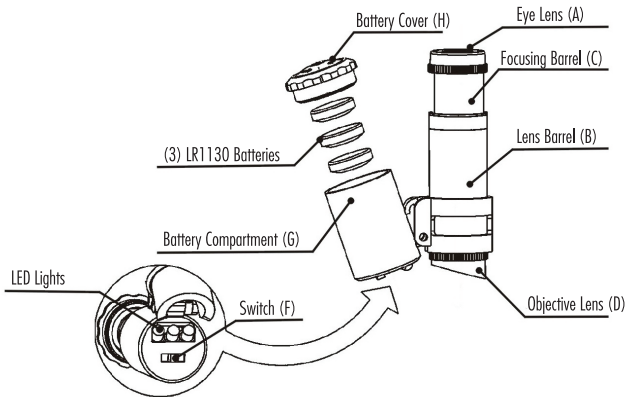

BATTERY REPLACEMENT:

1. Loosen the Battery Cover (H) by turning it counterclockwise.
2. Remove the old batteries from the Battery Compartment (G) and insert the new batteries.

Please Note: This microscope uses (3) LR1130 batteries.

3. Replace the Battery Cover (H) and tighten it by turning it clockwise.

DISTRIBUTED BY



DL WHOLESALE
DLWHOLESALE.COM