





LED SLIM-SCOPE 100X

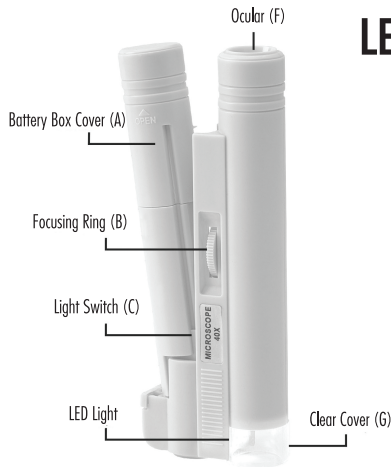

INSTRUCTIONS:

1. Open the Battery Box Cover (A) and insert (2) AA batteries (not included) with the positive ends pointing down, then close the Battery Box Cover (see diagram at bottom).
2. Place the sample to be viewed on a flat surface.
3. Place the microscope on the sample with the Clear Cover (G) facing downwards, towards the sample.
4. If needed, turn on the LED light using the Light Switch (C).
5. Look through the Ocular (F) and use the Focusing Ring (B) to adjust focus and clarity.

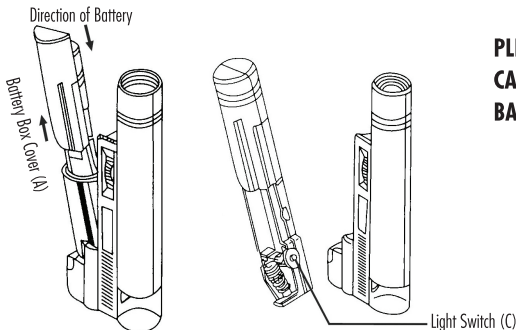

PLEASE NOTE: THIS MICROSCOPE REQUIRES (2) AA CARBON BATTERIES. PLEASE DO NOT USE ALKALINE BATTERIES, AS THEY MAY NOT FIT.

DISTRIBUTED BY



DL WHOLESALE
DLWHOLESALE.COM